

Regarding the use of clamps on the YZ1515X 6 rollers pump head, see the following instructions for details.

Installation method of the clamps

1. After installing the tube to the proper position of the pump head, open the clamp by hand gently, and clamp it to a position about 1cm on both sides of the pump head on the tube.
2. Use your hands or needle-nose pliers to clamp the clamp to the corresponding number of teeth in the instructions.



Not clamped



Clamping two teeth

3. Push the clamps to the state close to the pump head.


The clamp model, usage quantity (shipping quantity=single-side usage quantity×2) and number of clamping teeth corresponding to each type of tube are shown in the following table:

1. Tube Clamp Numbers and Clamping Teeth Number

Silicone Tube	Tube holder/Clamp	Numbers (single side)	Clamping teeth number
13#	Tube holder and clamp	2 holders and 1 clamp	1
14#	14# (3.8-4.3mm) Clamp	1	5
19#	19# (4.4-4.9mm) Clamp	1	1
16#	16# (5.0-5.5mm) Clamp	1	4
25#	25# (6.1-6.5mm) Clamp	2	3
17#	17# (7.6-8.5mm) Clamp	2	4
18#	18# (9.1-10.2mm) Clamp	2	4

Pharmed Tube	Tube holder/Clamp	Numbers(single side)	Clamping Teeth number
13#	Tube holder and clamp	1 holder and 1 clamp	1
14#	14# (3.8-4.3mm) Clamp	1	5
19#	19# (4.4-4.9mm) Clamp	1	1
16#	16# (5.0-5.5mm) Clamp	1	3
25#	25# (6.1-6.5mm) Clamp	1	1
17#	17# (7.6-8.5mm) Clamp	1	3
18#	18# (9.1-10.2mm) Clamp	2	4

Note: The number of teeth mentioned in the table above refers to the number of teeth on the outside of the clamp.

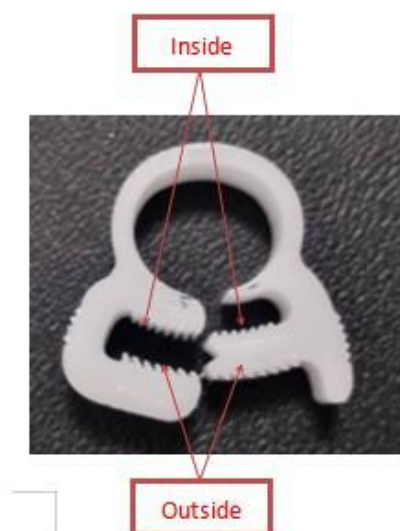

2. Picture for 13# tube with holder and clamp

